

## CE-MHC Class I Control Peptide Pool

**Catalog #:** CTL-CE-003, PA-CEF-003

**Lot:** \_\_\_\_\_

**Product:** CE-MHC Class I Control Peptide Pool

**Formulation:** Each vial contains a total of 400µg lyophilized peptide powder consisting of 20 individual peptides, 20µg of each peptide.

**Description:** The CE-MHC Control Peptide Pool contains 20 peptides, each corresponding to a defined HLA class I-restricted T cell epitope from cytomegalovirus and Epstein-Barr virus.

**Activity:** The CE-Control Peptide Pool stimulates corresponding peptide-specific CD8+ memory T cells to release IFN-γ. These peptides are restricted by HLA -A2, -A3, -A11, -A24, -B7, -B8, -B18, -B27, -B35, and -B44 molecules, and have been shown to elicit recall responses in the majority of individuals expressing these rather common MHC class I alleles – most humans have been previously exposed to these pathogens. Because of promiscuous binding motifs, these peptides can occasionally bind other MHC class I, but not class II molecules, further extending the range of individuals in whom they elicit a CD8 cell recall response. Therefore, the majority of randomly selected human donors will respond to the CE-Control peptide Pool.


**Recommended Use:** The CE-Control Peptide Pool is recommended as the positive control for detecting antigen-specific CD8+ cells in human PBMC, for example, when performing cytokine assays for immune monitoring purposes. Such assays include IFN-γ measurements by ELISPOT and intracytoplasmic cytokine staining (ICS) for accurate frequency measurements of the cytokine producing CD8 cells, or cytokine bead arrays (CBA), and ELISA, for a semi-quantitative read out.

**Instruction for Use:** **Stock solution:** Flick tube to ensure all powder is at the bottom of the tube. Add 10µl tissue culture grade DMSO followed by 40µl of sterile double-distilled water. Vortex briefly and maintain sterile handling! Add 450µl of CTL-Test™ Medium (CTLT-005) and vortex briefly. Flick tube to ensure that all liquid is at the bottom of the tube. The stock solution (500µl, 20x) is ready for use. The stock can be stored at 4°C for one week. CTL recommends to aliquot and store the stock solution at -20° to -80°C for long-term storage.

CTL recommends using Peptide Pool at 2µg/ml of final peptide concentration, that is at 1:20 dilution of the stock. For example, prepare a 1:10 dilution of the stock in CTL-Test™ Medium (CTLT-005) as the working solution, and add it 1+1 (v/v) to the medium containing the PBMC. For ELISPOT assays, we recommend plating 100µl of this 2x peptide working solution per well, directly into the ELISPOT plate, followed by adding the 100µl of the PBMC for a 24h assay duration. For exact frequency measurements, we recommend testing in triplicate. Because serum is the greatest variable in assay performance, we recommend the use of serum-free media at all steps of the



## Product Data Sheet

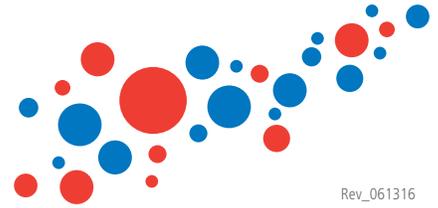

Rev\_061316

assay. CTL offers such Serum-free Media (CTL-Test™ Medium, CTLT-005) that has been customized for low-background/high-signal performance with PBMC. Investigators are advised to determine optimal cell concentration for individual applications for ELISPOT and related cytokine assays that is between 100,000 and 800,000 PBMC per well. CTL recommends 300,000 per well. MHC class I restricted peptides induce maximal CD8 cell stimulation when added 1+1 (vol+vol) to the PBMC.

**Storage:** In powder form at 4°C. The stock solution can be stored at 4°C for one week, however aliquoted and stored at -20° to -80°C is more suitable for storing long-term.

**Peptides:** CLGGLTMV (EBV, HLA-A2), GLCTLVAML (EBV, HLA-A2), NLVPMVATV (HCMV, HLA-A2), RVRAY-TYSK (EBV, HLA-A3), RLRAEAQVK (EBV, HLA-A3), AVFDRKSDAK (EBV, HLA-A11), IVTDFSVIK (EBV, HLA-A11), ATIGTAMYK (EBV, HLA-A11), DYCYNLNKEF (EBV, HLA-A24), RPPFIRRL (EBV, HLA-B7), RAKFKQLL (EBV, HLA-B8), FLRGRAYGL (EBV, HLA-B8), QAKWRLQTL (EBV, HLA-B8), SDEEEAIVAYTL (HCMV, HLA-B18), RRIYDLIEL (EBV, HLA-B27), YPLHEQHGM (EBV, HLA-B35), IPSINVHHY (HCMV, HLA-B35), EENLLDFVRF (EBV, HLA-B44), EFFWDANDIY (HCMV, HLA-B44), TPRVTGGGAM (HCMV, HLA-B7).

**References:** Currier JR, Kuta EG, Turk E, Earhart LB, Loomis-Price L, Janetzki S, Ferrari G, Birx DL, Cox JH. A panel of MHC class I restricted viral peptides for use as a quality control for vaccine trial ELISPOT assays. *J Immunol Methods* 260:157-172, 2002.

Mwau M, Michael A and Hanke T. Design and validation of an enzyme-linked immunospot assay for use in clinical trials of candidate HIV vaccines. *AIDS Res Hum Retroviruses* 18: 611-8, 2002.

**For laboratory research use only. Not for use in diagnostic or therapeutic procedures.**